

BIOTAQ™ DNA Polymerase

Shipping: On Dry/Blue Ice Catalog numbers
Batch No.: See vial BIO-21040: 500 Units (1 x 100 µL)
Concentration: 5 u/µL BIO-21060: 2500 Units (5 x 100 µL)



Store at -20°C

Storage and stability:

The BIOTAQ is shipped on dry/blue ice. On arrival store at -20 °C for optimum stability. Repeated freeze/thaw cycles should be avoided.

Expiry:

When stored under the recommended conditions and handled correctly, full activity of the kit is retained until the expiry date on the outer box label.

Safety precautions:

Please refer to the material safety data sheet for further information.

Unit definition:

One unit is defined as the amount of enzyme that incorporates 10nmoles of dNTPs into acid-insoluble form in 30 minutes at 72 °C.

Notes:

Research use only.

BIOTAQ, BioMix and HyperLadder are Trademarks of Bioline.

Features

- Premium Taq polymerase suited to a wide range of applications
- Amplifies fragments ≤5 kb
- Available as ready-to-use 2x reaction mixes (BioMix™/BioMix Red)

Applications

- Routine PCR applications
- TA cloning

Description

BIOTAQ™ is widely used by molecular biologists that have come to depend upon the robust performance of this reagent.

BIOTAQ is a highly purified thermostable DNA polymerase offering very high yield over a wide range of PCR templates, and is the ideal choice for most assays. BIOTAQ is a robust preparation and consistently delivers high yields with minimal background. BIOTAQ possesses 5'-3' exonuclease activity and leaves an 'A' overhang such that the PCR product is suitable for effective integration into TA cloning vectors.

BIOTAQ is supplied with 10x NH₄-based reaction buffer, which provides optimal conditions for most experiments. Additional MgCl₂ is provided to allow reaction conditions to be adjusted to suit the template.

Components:

Reagent	500 Units	2500 Units
BIOTAQ DNA Polymerase	1 x 100 µL	5 x 100 µL
10x NH ₄ Reaction Buffer	2 x 1.2 mL	10 x 1.2 mL
50mM MgCl ₂ Solution	1 x 1.2 mL	5 x 1.2 mL

General Considerations:

The optimum concentration of Mg²⁺ is 3 mM and should only be increased above this if absolutely necessary. For first tests, use no less than 2.5 units of BIOTAQ in a 50 µL reaction.

Citations: (http://www.bioline.com/h_scholar.asp)

1. Steinfurth, A., *et al. Endangered Species Res.* **39**: 293-302 (2019).
2. Murube, E., *et al. Crop and Pasture Sci.* **68(11)**:1006-1012 (2017).
3. Bennett, K.L., *et al. Mol. Ecol.* **25(17)**: 4337-4354 (2016).
4. Lutes, A.A. *et al. PNAS* **108**: 9910 - 9915 (2011).
5. Tucker, B.A. *et al. PNAS* **108**: E569 - E576 (2011)
6. Amaral, I.P. & Johnston, I.A. *J. Exp. Biol.* **214**: 2125-2139 (2011)
7. Coutinho, C.P., *et al. Infect. Immun.* **79**: 2950-2960 (2011)
8. Lora, J., *et al. PNAS* **108**: 5461-5465 (2011)
9. Levenberg, S. *et al. Nature Prot.* **5**: 1115-1126 (2010)
10. Tokuriki, N. & Tawfik, D.S. *Nature* **459**: 668-673 (2009)
11. Takada, S. & Mano, H. *Nature Prot.* **2**: 3136-3145 (2007)
12. López-Lluch, G., *et al. PNAS* **103(6)**: 1768-73 (2006)

PCR Reaction Conditions (for a 50 µL reaction)

10x NH ₄ Reaction Buffer	5 µL
50 mM MgCl ₂ Solution	3 µL
100 mM dNTP Mix (see below)	0.5 µL
Template and primers	As required
BIOTAQ	1 µL
Water (ddH ₂ O)	Up to 50 µL

Bioline 100 mM dNTP Mix is available as a separate product (Cat. No: BIO-39028)

Denature: 94-96 °C;
Extension: 70-72 °C allowing 15-30 seconds per kb

This data is intended for use as a guide only; conditions will vary from reaction to reaction and may need optimization.

Associated Products:

Product	Pack size	Cat. No.
dNTP Set	4 x 25 µmol	BIO-39025
dNTP Mix	500 µL	BIO-39028
HyperLadder™ 1kb	200 Lanes	BIO-33025

Bioline Reagents Ltd
UNITED KINGDOM

Tel: +44 (0)20 8830 5300
Fax: +44 (0)20 8452 2822

Bioline USA Inc.
USA

Tel: +1 901 382 8716
Fax: +1 901 382 0027

Bioline GmbH
GERMANY

Tel: +49 (0)3371 60222 00
Fax: +49(0)3371 60222 01

Bioline (Aust) Pty. Ltd
AUSTRALIA

Tel: +61 (0)2 9209 4180
Fax: +61 (0)2 9209 4763