Tube Roller

Instruction Manual (v.1)

N2400-7010

Tube Roller

CE

www.starlabgroup.com

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STARLAB certifies that

N2400-7010 Tube Roller

has been manufactured and inspected in accordance

with the following European standards:

2006/95/EC Low Voltage Directive

2004/108/EC Electromagnetic Compatibility Directive

Harmonized Standards:

EN 61010-1:2010

EN 61326-1:2006

As an ISO 9001 and ISO 13485 certified company,
our working procedures and processes are checked and audited regularly
and the quality of our products is continuously monitored.

STARLAB International GmbH

TUBE ROLLER

Cat. No. N2400-7010

1. Safety instructions

Before operation

- Read the Instruction Manual carefully before use.
- Ensure that only trained staff use the instrument. The instrument may only be opened by a qualified and trained technician.
- Make sure the socket is earthed before use.
- Set the instrument on a stable, clean, dry, non-slip and fireproof surface.
- Do not operate in hazardous areas.
- Before each use, check the instrument for damage. Do not use damaged components.
- The operation voltage noted on the label must match.
- Always disconnect the plug before fitting accessories.
- When working, wear personal protection to avoid the risk from splashing of liquids and possible glass breakage which can result from mechanical vibrations.

During operation

- Do not touch any moving parts during operation.
- If the instrument does not run smoothly decrease the motor speed.
- Set the "speed-control knob" to the lowest position first, then gradually increase the speed.

2. Product description

The instrument is designed for mixing liquids.

Normal use: The instrument is designed for use in professional laboratories. Do not use it in residential areas.

3. Inspection

Unpack the equipment carefully and check for any damages. If the instrument is faulty you should contact STARLAB or your STARLAB distributor promptly.

The packing includes:

- 1 Tube Roller N2400-7010
- 1 EU-power supply and 1 UK-power supply
- 1 Instruction manual

Please check all items are present.

4. Ambient conditions

Power supply: 100-240 V 50/60 Hz

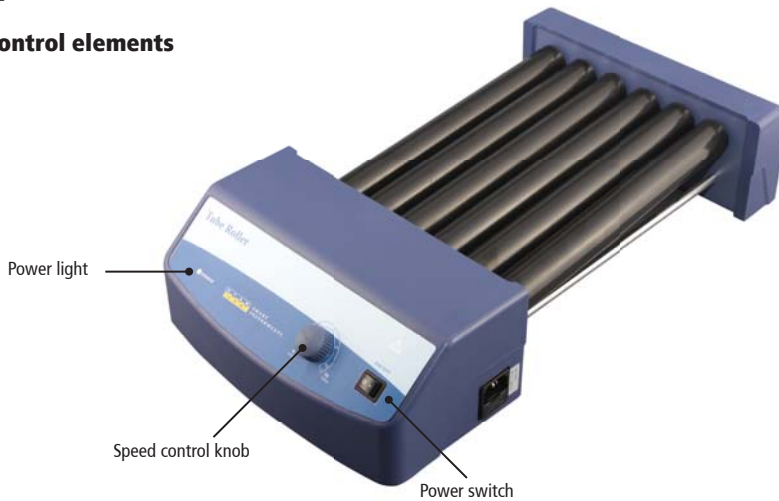
Ambient temperature: 4 – 40° C

Relative humidity: < 80 %

Place the instrument on a safe and stable surface.

5. Operation

5.1 Control elements

5.2 Operating

Switch on the instrument with the power switch. Turn the speed control knob to set the rated speed. Tubes can now be placed on the rolls. Switch off the instrument with the power switch and turn the speed control knob to "0".

6. Maintenance and cleaning

- Proper maintenance and cleaning will keep the instrument in good working order and will lengthen its life time.
- Be careful not spray cleaning fluid into the instrument. Unplug the power cord when cleaning.
- Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents the user must ascertain with the manufacturer that this will not destroy the instrument.
- Wear protective gloves during cleaning of the instrument.
- When not in use for long periods, switch off the instrument and store it in a dry, clean and stable place at room temperature.

Warranty

This instrument is under warranty to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty claim expires if the instrument is opened by untrained or unqualified staff. Our "General terms and conditions" apply.

Service

If help is needed, please contact STARLAB or your STARLAB distributor.
See www.starlabgroup.com for your local contact.

Please provide the customer care representative with the following information:

- Serial Number (on the rear panel)
- Description of problem (i.e., hardware or software)
- Methods and procedures adopted to resolve the problems
- Your contact information

7. Technical specifications

Parameter	Data
Voltage	100 – 240 V
Frequency	50/60 Hz
Power	25 W
Number of rollers	6
Size of rollers	280 x 30 mm
Max. Capacity	4 kg
Motor type	DC motor
Protection class	IP21
Weight	4,5 kg
Permitted ambient temperature	4 – 40 °C
Permitted relative humidity	< 80 %
Dimensions (WxDxH)	260 x 460 x 120 mm

8. Ordering information

Cat. No.	Description
N2400-7010	Tube Roller

Please visit our website www.starlabgroup.com, for details on our other instruments and our comprehensive range of consumables for the laboratory.

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STARLAB GmbH
STARLAB (UK), Ltd
STARLAB S.r.l.